

## Software Download and Installation

**Step 1:** Data logger software is available online at

<http://logtagrecorders.com/software/lta3/download/>

Ensure the LogTag cradle is **UNPLUGGED** before installing or updating software. **DO NOT** plug the cradle in until the software fully installs or updates. Failure to do this may result in disabled USB port.



**Step 2:** Click “Download Now”

**Step 3:** When the download is complete, click “open file” or “run”. If you are unable to find the file, look in your recent downloads.

**Step 4:** Agree to all “Terms & Conditions.” **Note:** You may need administrative privileges to download software onto your organization's computer.

## Software Settings (Configuration Part 1)


**Step 1:** Click on the icon on your desktop called LogTag Analyzer and open the program.

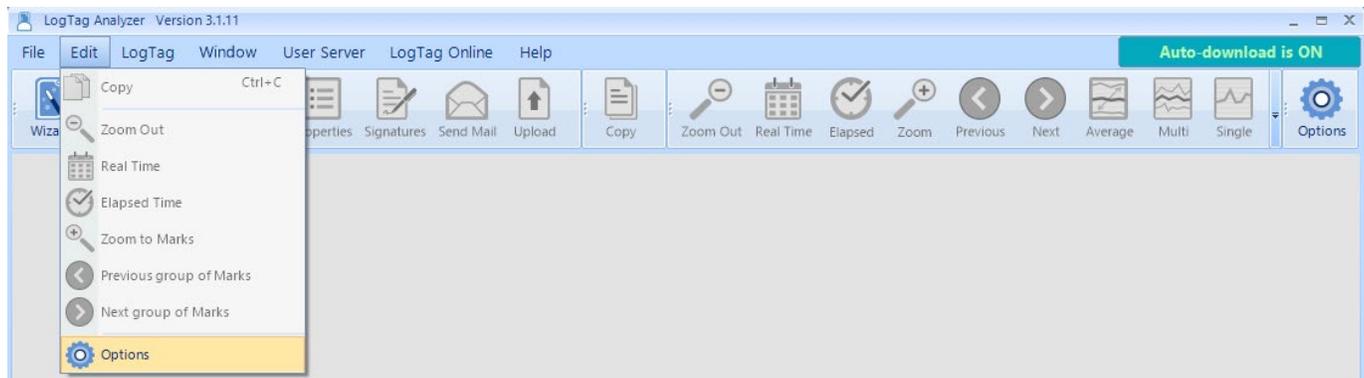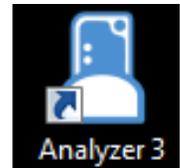


**Tip:** Some of the most common settings issues occur with reconfiguration, and location of the data file. See Automation and File and Folder Settings on page 4.

**Step 3:** Go through each of the below screens and update the software settings accordingly. **Do not click the “OK” button until all steps are complete.**


a.) **General Settings:** All temperatures should be in “Celsius”, and display “All Readings”

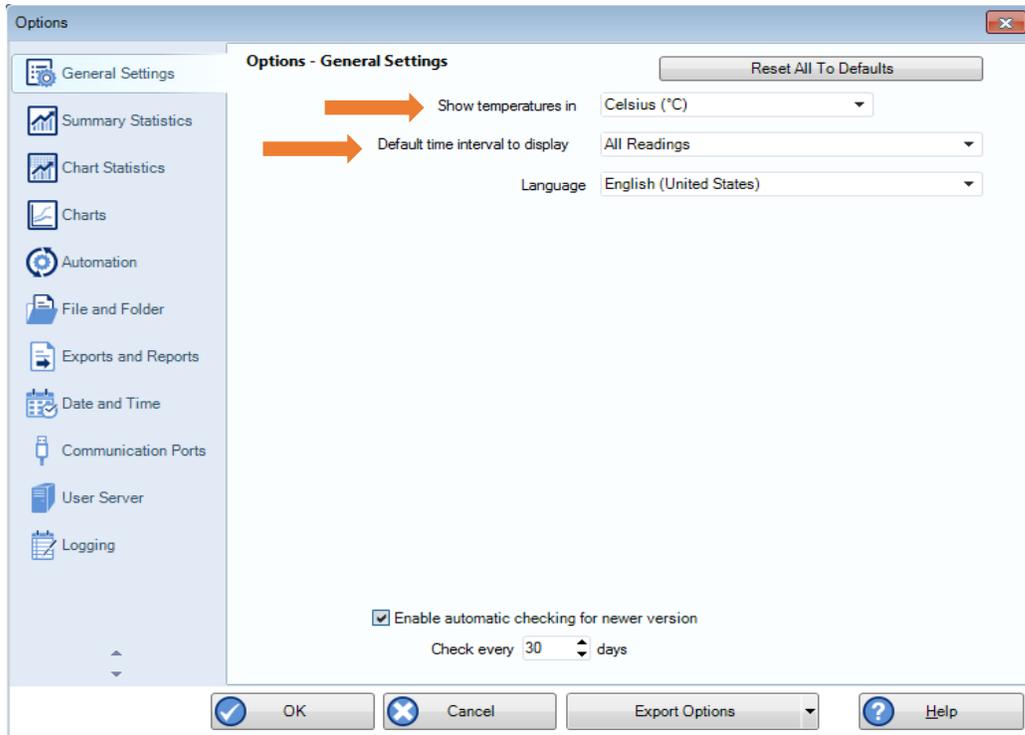


b.) **Summary Statistics:** Only the three boxes indicated should be checked off.

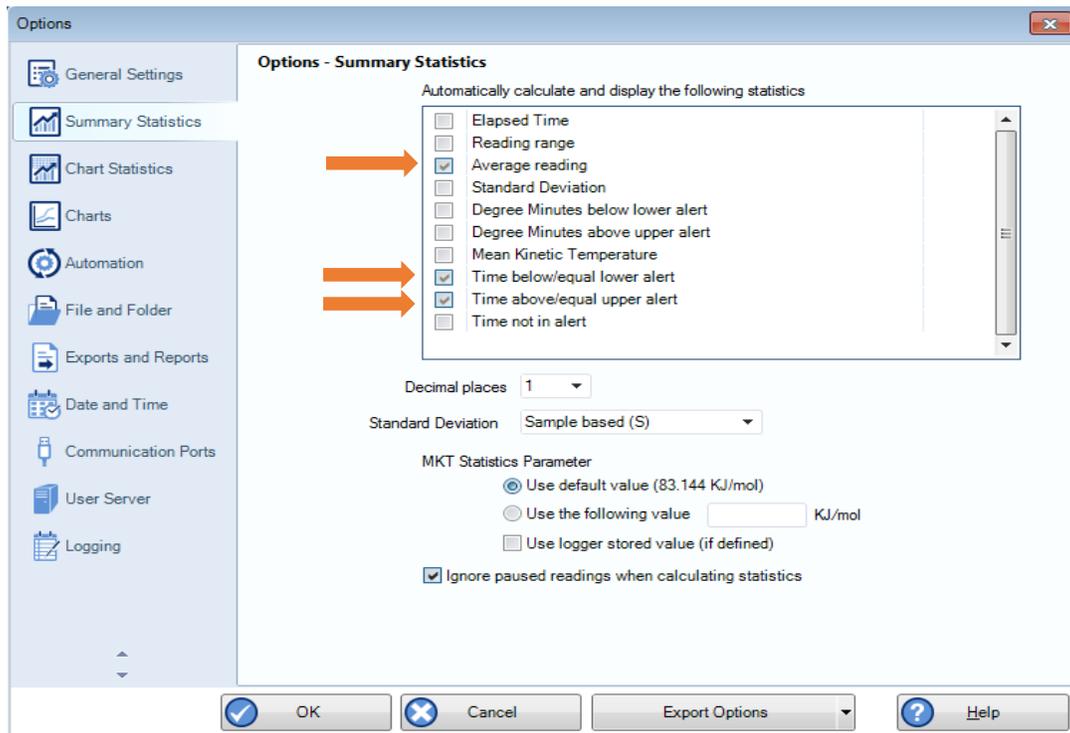


c.) **Chart Statistics:** Only the three boxes indicated should be checked off.

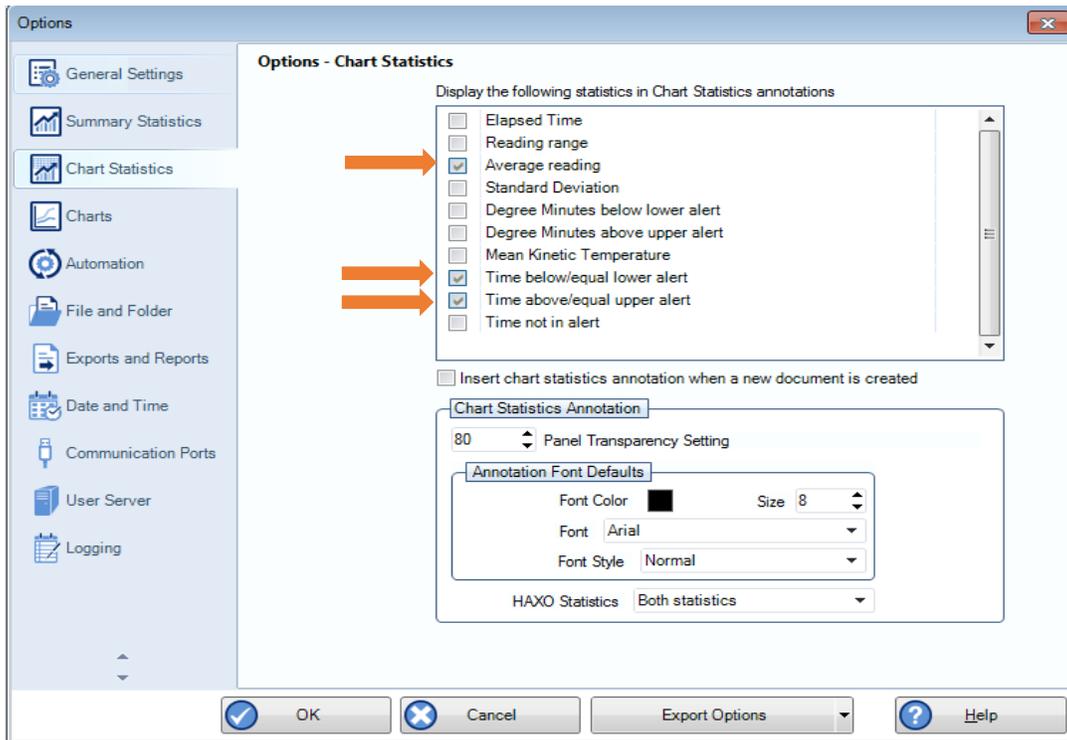

d.) **Charts:** Place a checkmark in each of the boxes indicated below. DO NOT change any of the other default settings.

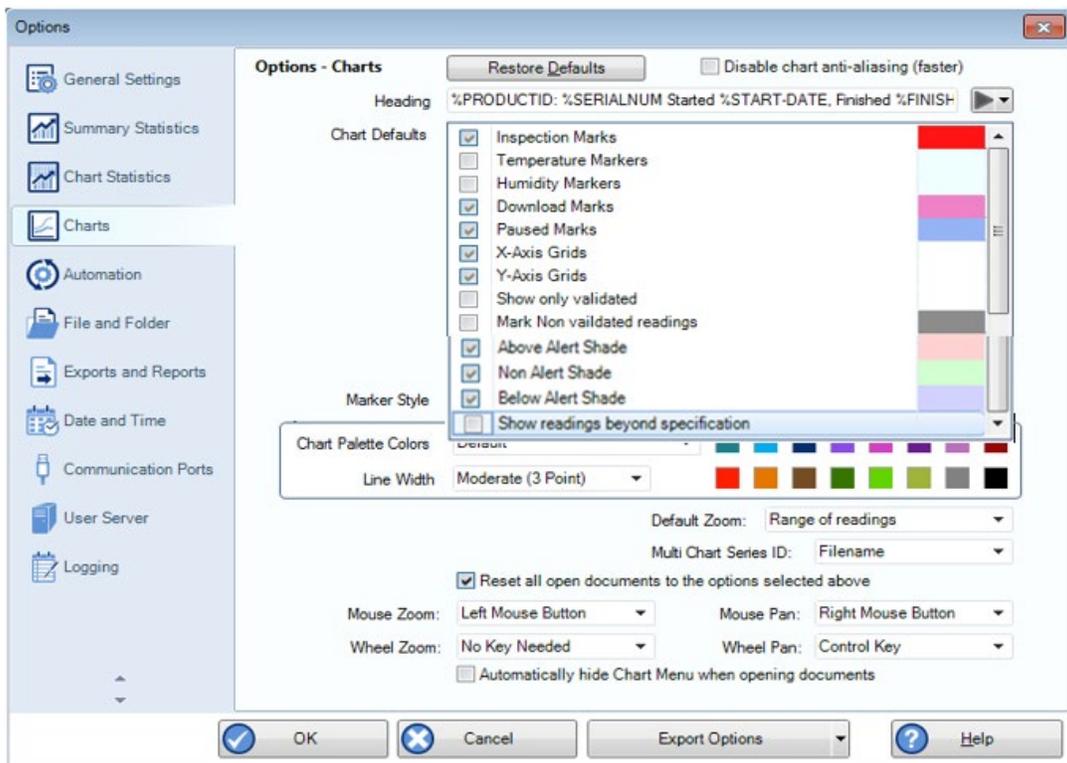


e.) Automation: **This is an important one.** Make sure the four boxes below are checked.

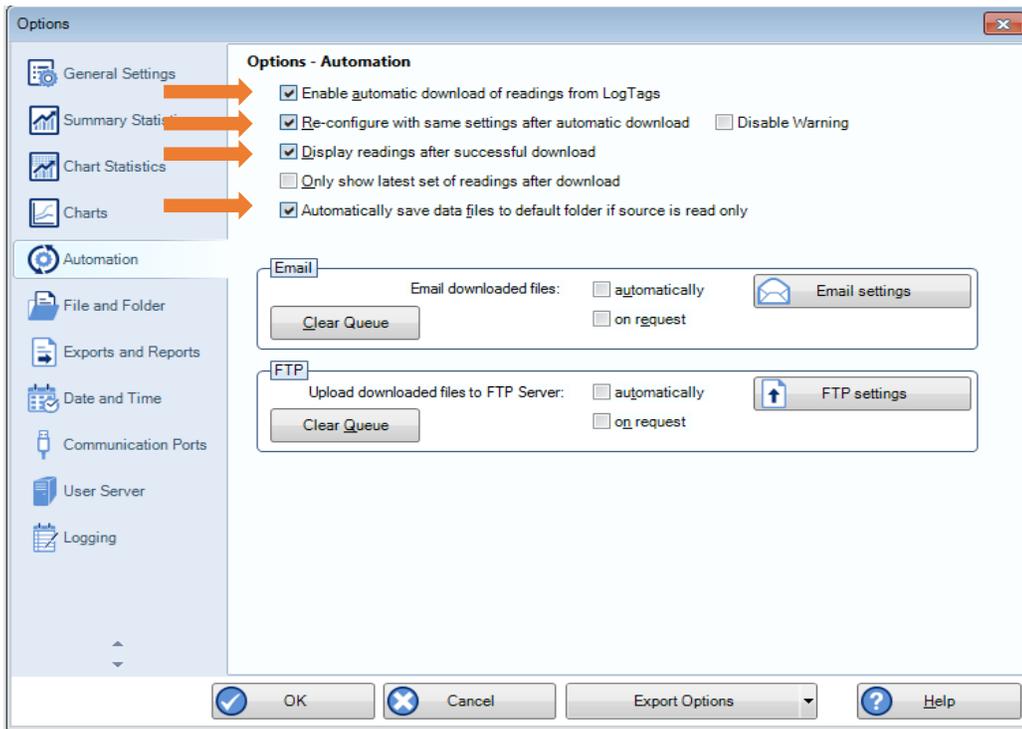


All four boxes must be checked for your logger to re-enter “READY” mode after download. If your settings are correct and your logger still does not reconfigure, please contact [AHS.VDHimmunizationProgram@vermont.gov](mailto:AHS.VDHimmunizationProgram@vermont.gov)

f.) File and Folder Settings: You do not need to change anything unless there is a specific folder the automatic downloads should go.

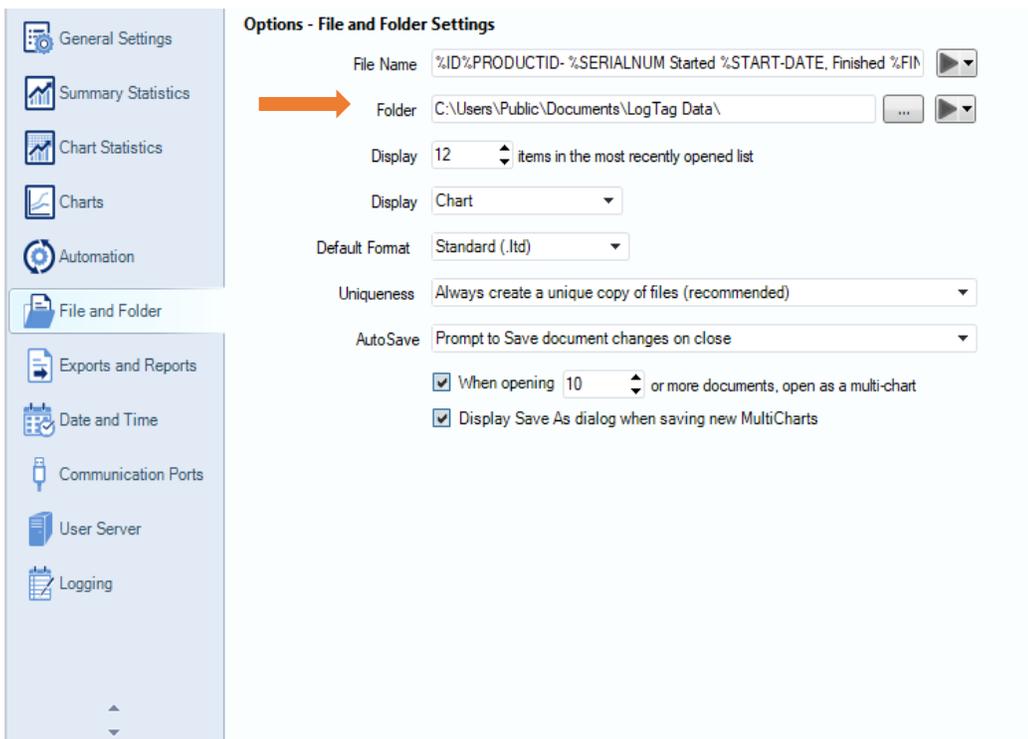


The software will default the save location to that shown. You may update this to whatever is convenient for your organization. If you have difficulty finding your data file, checking this setting may help.

g.) **Exports and Reports:** The only Export file type checked should be “LogTag File.”

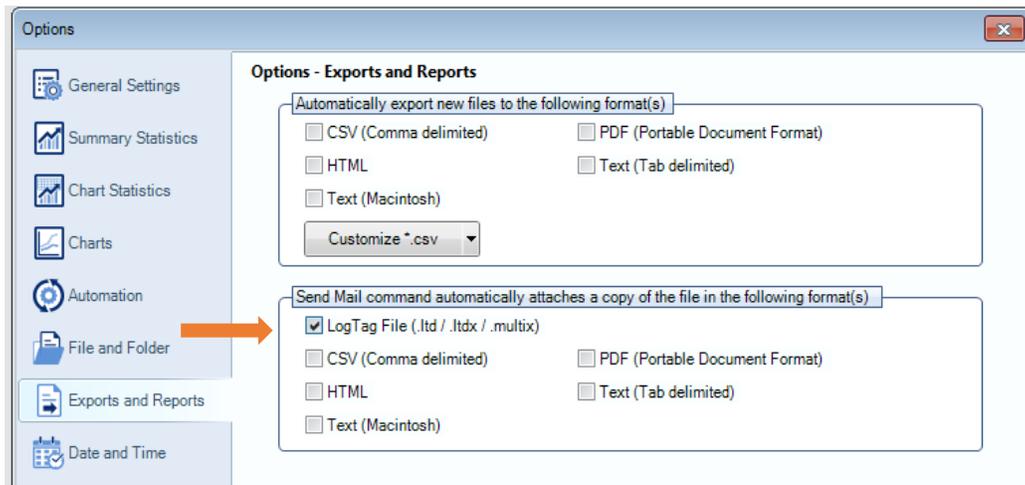

h.) **Date and Time:** No changes needed.

i.) **Communication Ports:** Check “USB Interface”

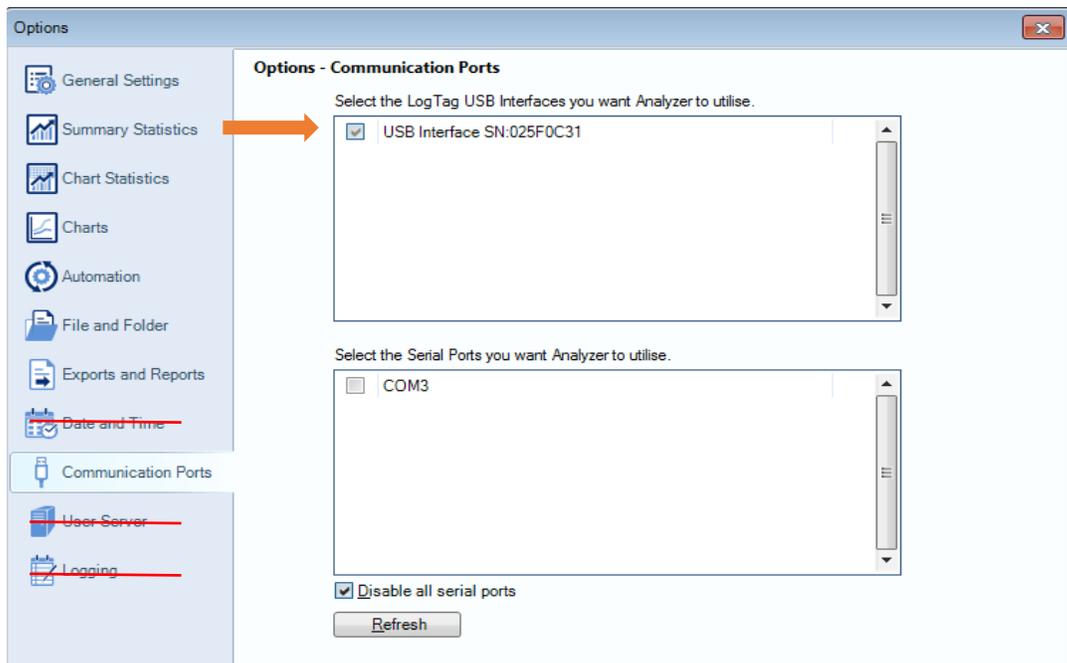

j.) **User Server:** No changes needed.

k.) **Logging:** No changes needed.

**Step 3:** Click “OK”

