

Reference Guide

Developed for
Vermont Department of Health



SensoScientific, Inc.

685 Cochran Street, Ste. #200, Simi Valley, CA 93065

www.sensoscientific.com

TABLE OF CONTENTS

- Cloud Portal Login 3
- Change Password 4
- Monitor Nodes 5
- Audit Nodes 6
- Confirm Alarms 7
- Turn Off Audible Alarm 8
- Audit Node Report 9
- Battery Status 10
- Contact Us 11


CLOUD PORTAL LOGIN

Use the SensoScientific Cloud Portal to view data logger records.

Initial Login

Log in to the
SensoScientific Cloud
by going to

cloud.sensoscientific.com

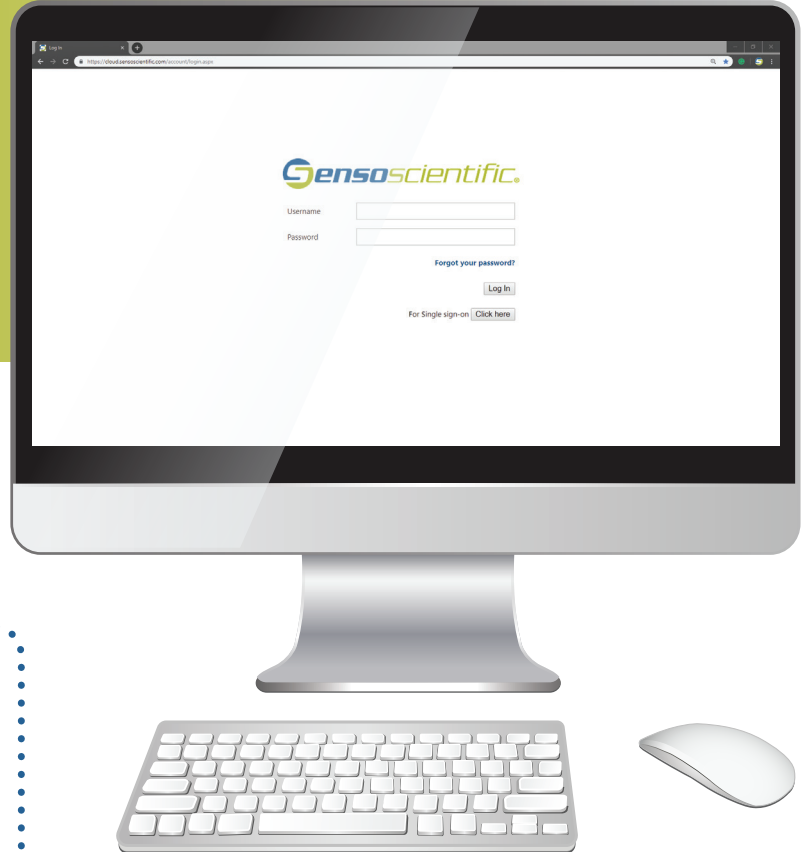

Please contact the **Vermont Immunization Program** for initial log in credentials and assistance. Keep in mind, *Passwords* and *Usernames* are **case-sensitive**.

CHANGE PASSWORD

Create new passwords.

- 1 Select the **“CONFIGURATION”** tab in the main menu.
- 2 Select the **“User”** tab in the submenu.
- 3 The submenu will reveal new options. Select the **“Change Password”** tab.
- 4 Fill out the fields and click the **“Change Password”** button.



1

ALARM | MONITORING | REPORT | **CONFIGURATION** | CALIBRATION | HELP

Company
Nodes
User
New User
Configure
Add/Remove Node
Change Password
Group
Settings
Notification

Change Password

2

You are about to change password for user : *demo*

Old Password :

New Password :

Confirm New Password :

3

Change Password

4

MONITOR NODES

View assigned nodes and check their current status.

- 1 Select the **“MONITORING”** tab in the main menu.
- 2 You will be automatically redirected to the **“Current Status”** page.
- 3 Each node will change color based on its **current status**.



ALARM | **MONITORING** | REPORT | CONFIGURATION | CALIBRATION | HELP

- Current Status
- Battery Status
- Reset Min/Max



Current Status



- GREEN:** All Values In Range
- YELLOW:** No Alarms - Value Out of Range, Not Yet Alarmed
- RED:** Unconfirmed Alarms Exist - Attention Required!
- PINK:** Suspended Alarm

Node Name	Node ID	Location	Alarm High	Alarm Low	Alarm Delay	Value	Graph	Last Updated	Value Status	Connectivity Status	Alarm Status	Notification Status
Lab Light	28329	Laboratory	N/A	N/A	60 (Mins)	No Light	24 Hrs	4/9/2019 2:23:42 PM	Out of Range	Connected	Alarmed	Active
99984E3DD	89949	Dietary	90 (°C)	-40 (°C)	30 (Mins)	83.7 (°C)	24 Hrs	4/9/2019 2:42:36 PM	In Range	Connected	Alarmed	Active
C8FD197B_1	81296	Room 5	90 (%RH)	-40 (%RH)	15 (Mins)	30 (%RH)	24 Hrs	4/9/2019 2:33:25 PM	In Range	Disconnected	Alarmed	Active
OR	11543	Pharmacy	8 (°C)	2 (°C)	0 (Mins)	7.6 (°C)	24 Hrs	4/9/2019 3:15:52 PM	In Range	Disconnected	Alarmed	Suspended
Temp 1	1245	Dietary	70 (%RH)	20 (%RH)	1 (Mins)	72 (%RH)	24 Hrs	4/9/2019 3:21:33 PM	Out of Range	Connected	Alarmed	Suspended
Temp 2	1247	Dietary	90 (°F)	-40 (°F)	30 (Mins)	21.1 (°F)	24 Hrs	4/9/2019 2:44:52 PM	In Range	Connected	No Alarms	Active
Walk In	86030	FoodService	46 (°F)	36 (°F)	30 (Mins)	30 (°F)	24 Hrs	4/9/2019 2:17:55 PM	In Range	Connected	No Alarms	Active
9884E3DD	1246	OR	85 (%RH)	30 (%RH)	60 (Mins)	60 (%RH)	24 Hrs	4/9/2019 2:21:28 PM	In Range	Connected	No Alarms	Active
Cafe	28329	NapaCafe	8 (°C)	2 (°C)	0 (Mins)	6 (°C)	24 Hrs	4/9/2019 3:44:42 PM	In Range	Connected	No Alarms	Active
Lab Ambient	81473	Laboratory	90 (°F)	-40 (°F)	30 (Mins)	21.1 (°F)	24 Hrs	4/9/2019 3:07:11 PM	In Range	Connected	No Alarms	Active
Vaccine	86027	Laboratory	30 (°C)	10 (°C)	90 (Mins)	28 (°C)	24 Hrs	4/9/2019 2:12:44 PM	Out of Range	Disconnected	No Alarms	Active
Visitor	86032		45 (°F)	35 (°F)	15 (Mins)	42.5 (°F)	24 Hrs	4/9/2019 2:15:46 PM	In Range	Disconnected	No Alarms	Active
Cafe	28329	NapaCafe	8 (°C)	2 (°C)	0 (Mins)	5.2 (°C)	24 Hrs	4/9/2019 3:17:57 PM	In Range	Connected	No Alarms	Active
Lab 2	81473	Laboratory	90 (°F)	-40 (°F)	30 (Mins)	21.1 (°F)	24 Hrs	4/9/2019 2:42:36 PM	In Range	Connected	No Alarms	Active

AUDIT NODES

Confirm that you have viewed a node's current status.

- 1 Select the **"MONITORING"** tab in the main menu.
- 2 Mark the checkbox in the rightmost column of the node(s) you want to audit.
- 3 Click the **"Audit Node"** button at the bottom of the page.



ALARM | **MONITORING** | REPORT | CONFIGURATION | CALIBRATION | HELP

- Current Status
- Battery Status
- Reset Min/Max

Current Status



Node Name	Node ID	Location	Alarm High	Alarm Low	Alarm Delay	Value	Graph	Last Updated	Value Status	Connectivity Status	Alarm Status	Notification Status	<input checked="" type="checkbox"/>
Lab Light	28329	Laboratory	N/A	N/A	60 (Mins)	No Light	24 Hrs	4/9/2019 2:23:42 PM	Out of Range	Connected	Alarmed	Active	<input checked="" type="checkbox"/>
99984E3DD	1249	Dietary	90 (°C)	-40 (°C)	30 (Mins)	83.7 (°C)	24 Hrs	4/9/2019 2:42:36 PM	In Range	Connected	Alarmed	Active	<input checked="" type="checkbox"/>
CBFD197B_T	1245	Room 5	90 (%RH)	-40 (%RH)	15 (Mins)	30 (%RH)	24 Hrs	4/9/2019 2:33:25 PM	In Range	Disconnected	Alarmed	Active	<input checked="" type="checkbox"/>
OR	11545	Pharmacy	8 (°C)	2 (°C)	0 (Mins)	7.6 (°C)	24 Hrs	4/9/2019 3:15:52 PM	In Range	Disconnected	Alarmed	Suspended	<input checked="" type="checkbox"/>
Temp 1	1247	Dietary	70 (%RH)	20 (%RH)	1 (Mins)	72 (%RH)	24 Hrs	4/9/2019 3:21:33 PM	Out of Range	Connected	Alarmed	Suspended	<input checked="" type="checkbox"/>
Temp 2	86025	Dietary	90 (°F)	-40 (°F)	30 (Mins)	21.1 (°F)	24 Hrs	4/9/2019 2:44:52 PM	In Range	Connected	No Alarms	Active	<input checked="" type="checkbox"/>
Walk In	46422	FoodService	46 (°F)	36 (°F)	30 (Mins)	30 (°F)	24 Hrs	4/9/2019 2:17:55 PM	In Range	Connected	No Alarms	Active	<input checked="" type="checkbox"/>
9884E3DD	86029	OR	85 (%RH)	30 (%RH)	60 (Mins)	60 (%RH)	24 Hrs	4/9/2019 2:21:28 PM	In Range	Connected	No Alarms	Active	<input checked="" type="checkbox"/>
Cafe	86030	NapaCafe	8 (°C)	2 (°C)	0 (Mins)	6 (°C)	24 Hrs	4/9/2019 3:44:42 PM	In Range	Connected	No Alarms	Active	<input checked="" type="checkbox"/>
Lab Ambient	89949	Laboratory	90 (°F)	-40 (°F)	30 (Mins)	21.1 (°F)	24 Hrs	4/9/2019 3:07:11 PM	In Range	Connected	No Alarms	Active	<input checked="" type="checkbox"/>
Vaccine	1246	Laboratory	30 (°C)	10 (°C)	90 (Mins)	28 (°C)	24 Hrs	4/9/2019 2:12:44 PM	In Range	Disconnected	No Alarms	Active	<input checked="" type="checkbox"/>
Visitor	1248		45 (°F)	35 (°F)	15 (Mins)	42.5 (°F)	24 Hrs	4/9/2019 2:15:46 PM	In Range	Disconnected	No Alarms	Active	<input checked="" type="checkbox"/>
PH Humid	11543	NapaCafe	8 (°C)	2 (°C)	30 (Mins)	5.2 (°C)	24 Hrs	4/9/2019 3:17:57 PM	In Range	Connected	No Alarms	Active	<input checked="" type="checkbox"/>
Lab 2	11544	Laboratory	90 (°F)	-40 (°F)	30 (Mins)	21.1 (°F)	24 Hrs	4/9/2019 2:42:36 PM	In Range	Connected	No Alarms	Active	<input checked="" type="checkbox"/>
Visitor 1	17931	Corporate	8 (°C)	2 (°C)	15 (Mins)	6 (°C)	24 Hrs	4/9/2019 3:44:42 PM	In Range	Connected	No Alarms	Active	<input checked="" type="checkbox"/>
Visitor 2	28327	Corporate	90 (°F)	-40 (°F)	30 (Mins)	21.1 (°F)	24 Hrs	4/9/2019 3:07:11 PM	In Range	Connected	No Alarms	Active	<input checked="" type="checkbox"/>
9984E3D_T	28328	NapaCafe	13.3 (°C)	12 (°C)	0 (Mins)	22.1 (°C)	24 Hrs	4/9/2019 2:22:56 PM	Out of Range	Disconnected	No Alarms	Active	<input checked="" type="checkbox"/>

Audit Node

CONFIRM ALARMS

Document the cause of alarms.

- 1 Select the “**MONITORING**” tab in the main menu.
- 2 Click the “**Alarmed**” button of the desired node. *(Not Displayed)*
- 3 Click the “**Take Action**” button to confirm the alarm. *(Not Displayed)*
- 4 Fill out the fields and click the “**Confirm**” button.

Sensoscientific.

1

ALARM | MONITORING | REPORT | CONFIGURATION | CALIBRATION | HELP

Diagnose Alarm

Diagnose Alarm

Node Name	Node ID	Node Type	Alarm Date	Alarm Type
Lab Light	28329	Light Sensor	4/8/2019 7:05:51 PM	Data Alarm

Root Cause:

Detail Description (optional)

Corrective Action (Max 500 characters)

Is the equipment working properly? Yes No

Confirm

4

If the vaccine storage unit is not working properly, you have the option to **suspend the node**. This will stop alarm notifications for the designated duration.

TURN OFF AUDIBLE ALARM

Deactivate a node's audible alarm.

- 1 Select the **"MONITORING"** tab in the main menu.
- 2 Click the **speaker icon** of the node you want to silence.



ALARM | **MONITORING** | REPORT | CONFIGURATION | CALIBRATION | HELP

Current Status

Battery Status

Reset Min/Max

Current Status



Audible Alarm: **Active**



Audible Alarm: **Inactive**



Node Name	Node ID	Location	Alarm High	Alarm Low	Alarm Delay	Value	Graph	Last Updated	Value Status	Connectivity Status	Alarm Status		<input type="checkbox"/>
Lab Light	28329	Laboratory	N/A	N/A	60 (Mins)	No Light		4/9/2019 2:23:42 PM	Out of Range	Connected	Alarmed		<input type="checkbox"/>
99984E3DD	1249	Dietary	90 (°C)	-40 (°C)	30 (Mins)	83.7 (°C)		4/9/2019 2:42:36 PM	In Range	Connected	Alarmed		<input type="checkbox"/>
C8FD197B_T	1245	Room 5	90 (%RH)	-40 (%RH)	15 (Mins)	30 (%RH)		4/9/2019 2:33:25 PM	In Range	Disconnected	Alarmed	N/A	<input type="checkbox"/>
OR	11545	Pharmacy	8 (°C)	2 (°C)	0 (Mins)	7.6 (°C)		4/9/2019 3:15:52 PM	In Range	Disconnected	Alarmed		<input type="checkbox"/>
Temp 1	1247	Dietary	70 (%RH)	20 (%RH)	1 (Mins)	72 (%RH)		4/9/2019 3:21:33 PM	Out of Range	Connected	Alarmed		<input type="checkbox"/>
Temp 2	86025	Dietary	90 (°F)	-40 (°F)	30 (Mins)	21.1 (°F)		4/9/2019 2:44:52 PM	In Range	Connected	No Alarms	N/A	<input type="checkbox"/>
Walk In	46422	FoodService	46 (°F)	36 (°F)	30 (Mins)	30 (°F)		4/9/2019 2:17:55 PM	In Range	Connected	No Alarms		<input type="checkbox"/>
9884E3DD	86029	OR	85 (%RH)	30 (%RH)	60 (Mins)	60 (%RH)		4/9/2019 2:21:28 PM	In Range	Connected	No Alarms	N/A	<input type="checkbox"/>
9884E3DD	86029	OR	85 (%RH)	30 (%RH)	60 (Mins)	60 (%RH)		4/9/2019 2:21:28 PM	In Range	Connected	No Alarms	N/A	<input type="checkbox"/>
9884E3DD	86029	OR	85 (%RH)	30 (%RH)	60 (Mins)	60 (%RH)		4/9/2019 2:21:28 PM	In Range	Connected	No Alarms		<input type="checkbox"/>
9884E3DD	86029	OR	85 (%RH)	30 (%RH)	60 (Mins)	60 (%RH)		4/9/2019 2:21:28 PM	In Range	Connected	No Alarms		<input type="checkbox"/>

AUDIT NODE REPORT

Generate a report of all audited nodes.

- 1 Select the **“REPORT”** tab in the main menu.
- 2 Select the **“Audit Node”** tab in the submenu.
- 3 Specify the parameters and click the **“Show Report”** button.



ALARM | MONITORING | **REPORT** | CONFIGURATION | CALIBRATION | HELP

Audit Node

- Graph Report
- Daily Report
- Weekly Report
- Advanced Report
- Corrective Action Report
- Performance Report
- Two Point Monitoring
- Mean Kinetic Temp
- Audit Node**
- Admin Audit Report
- Audit User Action
- Connection Report
- Low Battery Report
- Detail Graph Report
- Suspended Action



All Nodes Select Node
 Select Start Date Select End Date
Show Report



BATTERY STATUS

Check the battery status of your assigned nodes.

- 1 Select the “**MONITORING**” tab in the main menu.
- 2 Select the “**Battery Status**” tab in the submenu.
- 3 Each node will change color based on its current **battery status**.



ALARM | **MONITORING** | REPORT | CONFIGURATION | CALIBRATION | HELP

- Current Status
- Battery Status**
- Reset Min/Max



Battery Status



GREEN: High Battery - No Action Needed



YELLOW: Normal Battery - Extra Batteries Recommended



RED: Low Battery - Order Batteries

Node Name	Node ID	Location	Node Type	Status	Message	Last Updated	Alarm Status	Power Source
Lab Ambient	28327	Laboratory	Temperature		High	4/19/2019 11:38:48 AM	No Alarms	External Power
OR Pharmacy	46422	Pharmacy	Humidity		High	4/19/2019 11:34:28 AM	No Alarms	Internal Battery
Walk In Freezer	86025	Food Service	Temperature		High	6/26/2018 12:51:08 PM	No Alarms	N/A
Blood Bank	1245	Blood Bank	Temperature		High	4/19/2019 11:38:48 AM	No Alarms	External Power
Pharmacy Main	1247	Pharmacy	Temperature		Normal	4/19/2019 11:34:28 AM	No Alarms	Internal Battery
Specimen Fridge	11543	Laboratory	Humidity		Normal	6/26/2018 12:51:08 PM	No Alarms	N/A
Walk In Fridge 2	17931	Food Service	Temperature		High	4/19/2019 11:38:48 AM	No Alarms	External Power
Cafe Refrigerator	46423	NapaCafe	Temperature		High	4/19/2019 11:34:28 AM	No Alarms	Internal Battery
C8FD197B22EF_T	81296	Laboratory	Light		Low	6/26/2018 12:51:08 PM	No Alarms	N/A
Pharmacy Fridge	85236	Pharmacy	Temperature		High	4/19/2019 11:38:48 AM	No Alarms	External Power
Vaccine Sensor	28329	Corporate Office	Temperature		High	4/19/2019 11:34:28 AM	No Alarms	Internal Battery
Visitor Center	11544	NapaCafe	Temperature		High	6/26/2018 12:51:08 PM	No Alarms	N/A

CONTACT US

Vermont Department of Health Immunization Program

108 Cherry Street
Burlington, VT 05401

P. (802) 863-7638 or 1-800-640-4374
ahs.vdhimmunizationprogram@vermont.gov